

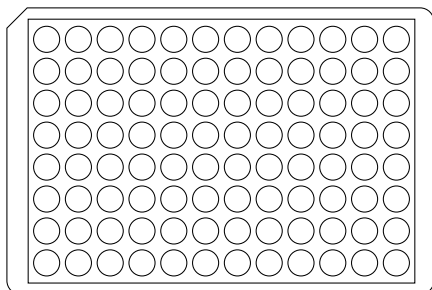


Magnetic Separator - 96 Well Plate 472236

Gather Components

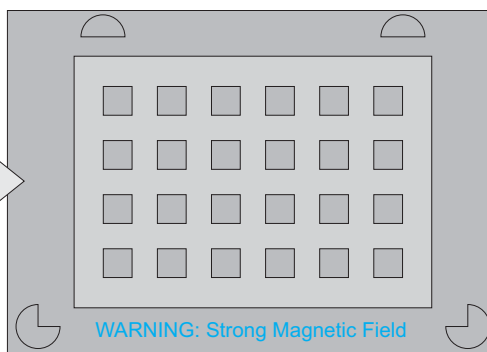


96-Well Plate
Each well containing a suspension of magnetic particles.

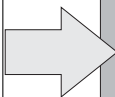

Magnetic Separator


1. Placement

Place the 96-Well Plate on to the Magnetic Separator.

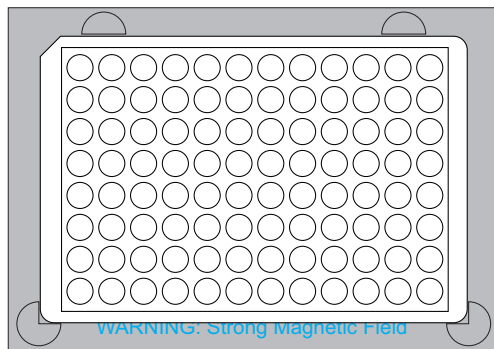

2. Empty the 96-Well Plate

While firmly holding the 96-Well Plate against the Magnetic Separator, invert over a sink to discard the liquid. The magnetic particles will be retained in the 96-Well Plate.

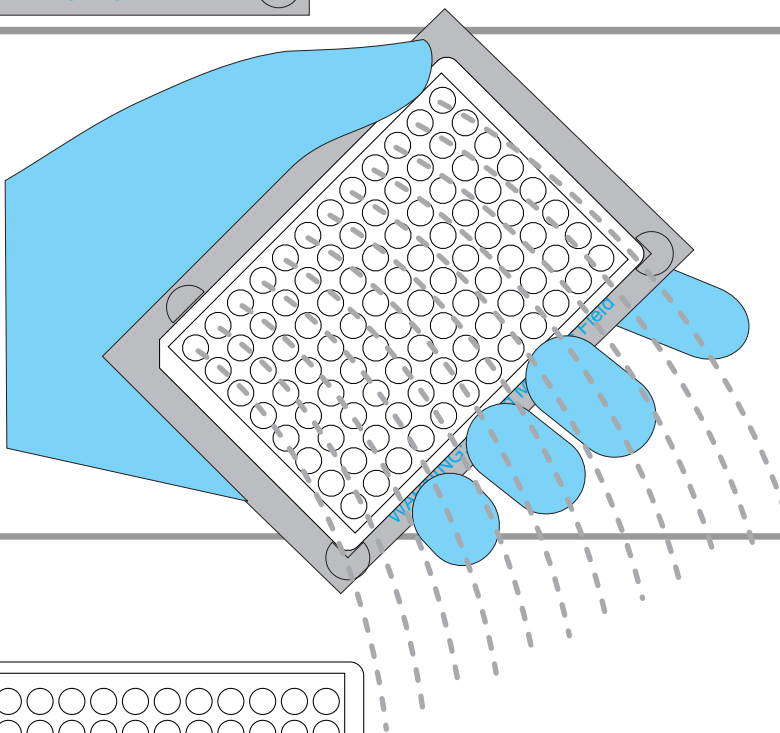

3. Remove the 96-Well Plate

The magnetic particles retained in the plate are now ready for further use.

