


Deoxynivalenol/Vomitoxin ELISA Plate 53016B


1. Addition of Enzyme Conjugate

Add 50 uL of the enzyme conjugate to the individual wells successively using a multi-channel pipette or a stepping pipette.

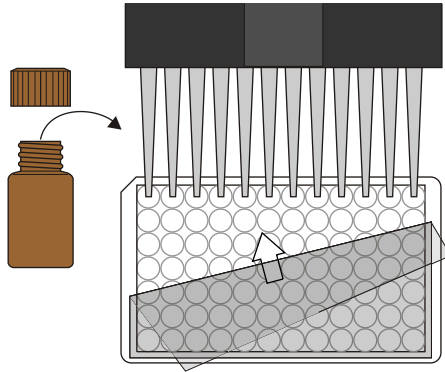


5. Addition of Substrate/Color Solution

Add 100 uL of substrate/color solution to the individual wells successively using a multi-channel pipette or a stepping pipette. cover the wells with parafilm or tape and mix the contents by moving the strip holder in a rapid circular motion on the benchtop. Be careful not to spill contents. Incubate the strips for 5 minutes at room temperature.

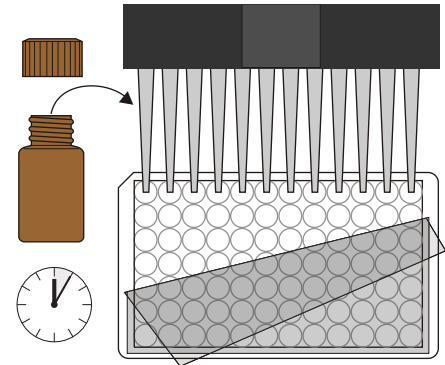


2. Addition of Standards, Samples

Add 50 uL of the standard solutions and the samples into the wells of the test strips. We recommend using duplicates or triplicates.

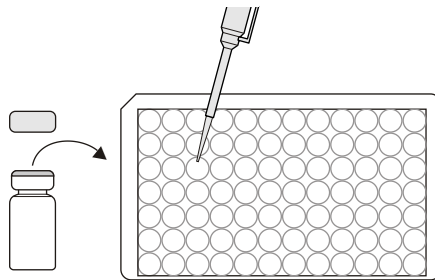


6. Addition of Stopping Solution

Add 100 uL of stop solution to the wells in the same sequence as for the substrate solution using a multi-channel pipette or a stepping pipette.

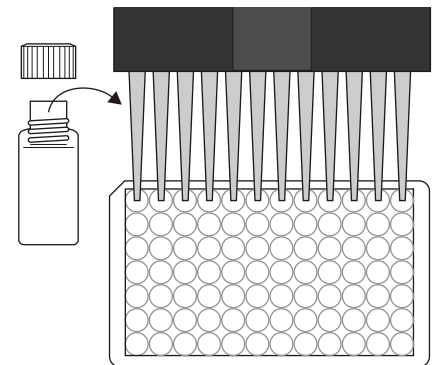


3. Addition of Antibody Solution

Add 50 uL of the antibody solution into the wells of the test strips. Then incubate at room temperature (15 - 30°C) for 10 minutes.

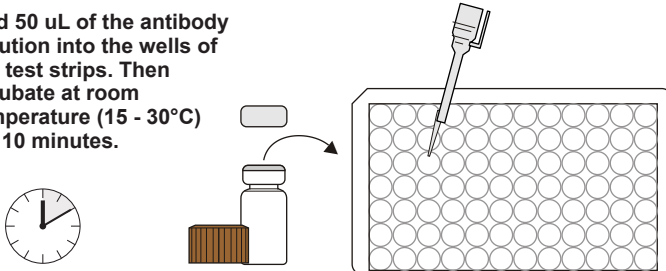

7. Measurement of Color

Read the absorbance at 450nm using a microplate ELISA reader. Calculate the results.

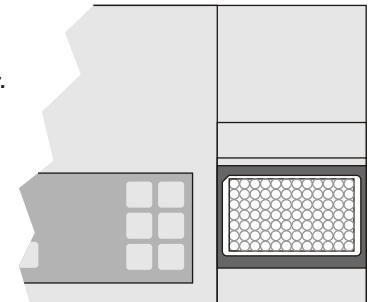

4. Washing of Plates

After incubation, remove the covering and vigorously shake the contents of the wells into a sink. Wash the strips five times with a multi-channel pipette or wash bottle using the 1X washing buffer solution. Please use at least a volume of 250 uL of washing buffer for each well and each washing step. Remaining buffer in the wells should be removed by patting the plate dry on a stack of paper towels.

